



< Example of DNA extraction >

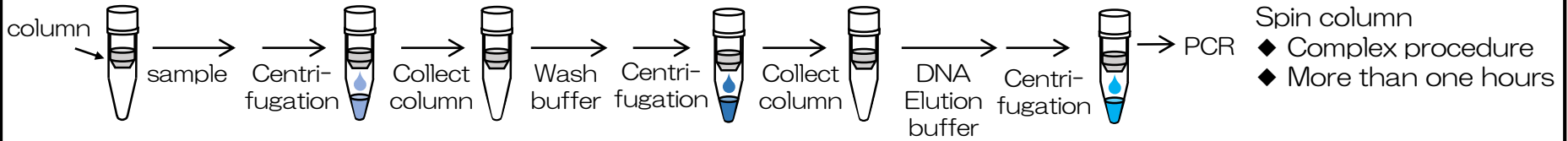

< Kaneka Easy DNA Extraction Kit version 2* >

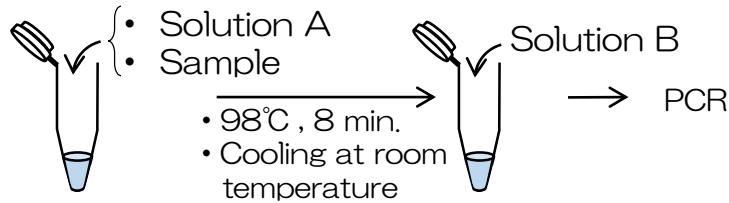

- Simple protocol
- 10 minute extraction

*This product is for research use only.

< Performance comparison with a spin column DNA purification kit >

Sample: bird feces.

Target sequence: mitochondria COI (648 bp)


<Performance comparison>

PCR following microchip electrophoresis

“Kaneka DNA Easy Extraction Kit version 2” can extracted DNA from bird feces at roughly the same efficiency level as a spin column kit.

Kaneka Easy DNA Extraction Kit version 2

PCR amplification (+) (+) (+) (-) (-) (+)

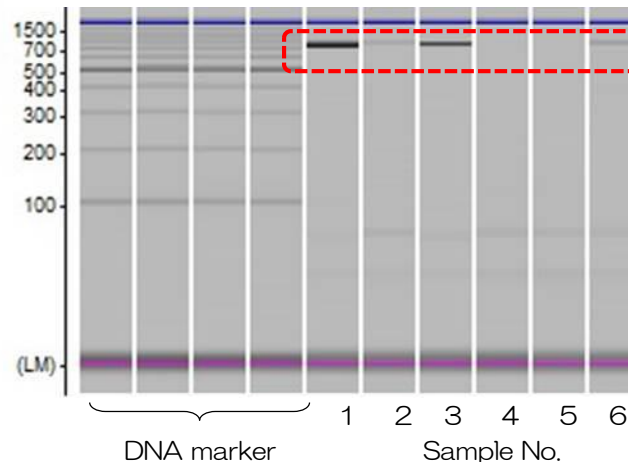

Spin column DNA purification kit (kit A)

PCR amplification (+) (+) (+) (-) (-) (+)

